

Laser Ablation Split Stream (LASS) Between Three Inductively Coupled Plasma Mass Spectrometers for Zircon Petrochronology

G. Craig¹, D. A. Frick², C. Bouman¹, N. Lloyd¹ and J. B. Schwieters¹

¹ Thermo Fisher Scientific (Bremen) GmbH, 28199 Bremen, Germany

² Helmholtz Centre Potsdam, GFZ German Research Centre for Geosciences, 14473 Potsdam, Germany

Overview

Purpose: To evaluate the effectiveness of laser ablation split stream (LASS) between three different inductively coupled plasma mass spectrometers (ICP-MS) for zircon geochronology.

Methods: LASS was achieved using disposable membrane pumps. Hf, U-Pb and rare earth element (REE) data were measured simultaneously by multicollector (MC), sector field (SF), and quadrupole (Q) ICP-MS respectively.

Results: Hf and U-Pb isotope ratios, and REE concentrations were determined successfully for four reference materials.

Introduction

In laser ablation analysis the choice of which type of ICP-MS to use is dependent on the application. Chronometers requiring epsilon level precision are measured on multicollector ICP-MS: for determination of the rare earth composition a quadrupole based instrument is sufficient. The development of LASS, where the laser aerosol is split between two ICP-MS, has allowed two isotopic (or trace elemental) systems to be measured simultaneously on a single ablation site. A primary LASS application is zircon and monazite petrochronology¹, for which two separate chronometers, U-Pb and Hf, and REE concentration, are often collected. By determining both the chronometers and the elemental composition for a single ablation site, a direct link between each can be established.

LASS analysis of common reference zircons and monazites combining three different Thermo Scientific™ ICP-MS, the Neptune Plus™ (for Hf), Element XR™ (for U-Pb) and iCAP™ RQ (for REE) is reported. The laser ablation gas stream was split between the three ICP-MS, following the approach was outlined by Frick *et al.*

Methods

Sample Reference Materials

- NIST SRM® 610
- 4 reference zircons and monazites
 - 91500
 - Plesovice
 - GJ-1
 - Mud Tank
- MUN artificial zircon - external Yb correction of Hf.

Laser Ablation

The LA system used was a Teledyne Photon Machines Analyte G2™ equipped with a HelEx II ablation cell.

- Spot Size – 50 μm
- Repetition Rate - 10 Hz
- Fluence – 4J cm⁻²
- Shot Count – 600
- He Outer Cell – 0.75 L min⁻¹
- He Cup Flow – 0.60 L min⁻¹

Thermo Scientific™ NEPTUNE Plus™

1.049 s integration time was used for all analyses. The Hf cup configurations was:

L4	L3	L2	L1	C	H1	H2	H3	H4
¹⁷¹ Yb	¹⁷³ Yb	¹⁷⁵ Lu	¹⁷⁶ Hf	¹⁷⁷ Hf	¹⁷⁸ Hf	¹⁷⁹ Hf		

Thermo Scientific™ Element XR™

Isotopes	²⁰² Hg	²⁰⁴ Pb	²⁰⁶ Pb	²⁰⁷ Pb	²⁰⁸ Pb	²³² Th	²³⁵ U	²³⁸ U
Segment Duration	5 ms	5 ms	26 ms	48 ms	26 ms	26 ms	26 ms	26 ms

Thermo Scientific™ iCAP™ RQ

Monitored Isotopes (Dwell time – 50 ms)

- ⁸⁹Y, ⁹¹Zr, ⁹³Nb, ¹³⁹La, ¹⁴⁰Ce, ¹⁴¹Pr, ¹⁴⁶Nd, ¹⁴⁷Sm, ¹⁵³Eu, ¹⁵⁷Gd, ¹⁵⁹Tb, ¹⁶³Dy, ¹⁶⁵Ho, ¹⁶⁹Tm, ¹⁷²Yb, ¹⁷⁵Lu, ¹⁷⁸Hf

LASS

Compact membrane pumps were used to perform each split, extracting ablated material from the central LA-MC-ICP-MS channel to the other ICP-MS. The membrane pumps extracted part of the ablation gas stream from a LA mixing chamber (Figure 1). The amount of material extracted could be varied for each pump.

- Split 1 – 0.75 L min⁻¹ pump at 88% efficiency.
- Split 2 – 0.40 L min⁻¹ pump at 94% efficiency.

Data Analysis

All data analysis was performed within Iolite v3.31³ (Igor Pro™ v6.37) using appropriate Data Reduction Schemes.

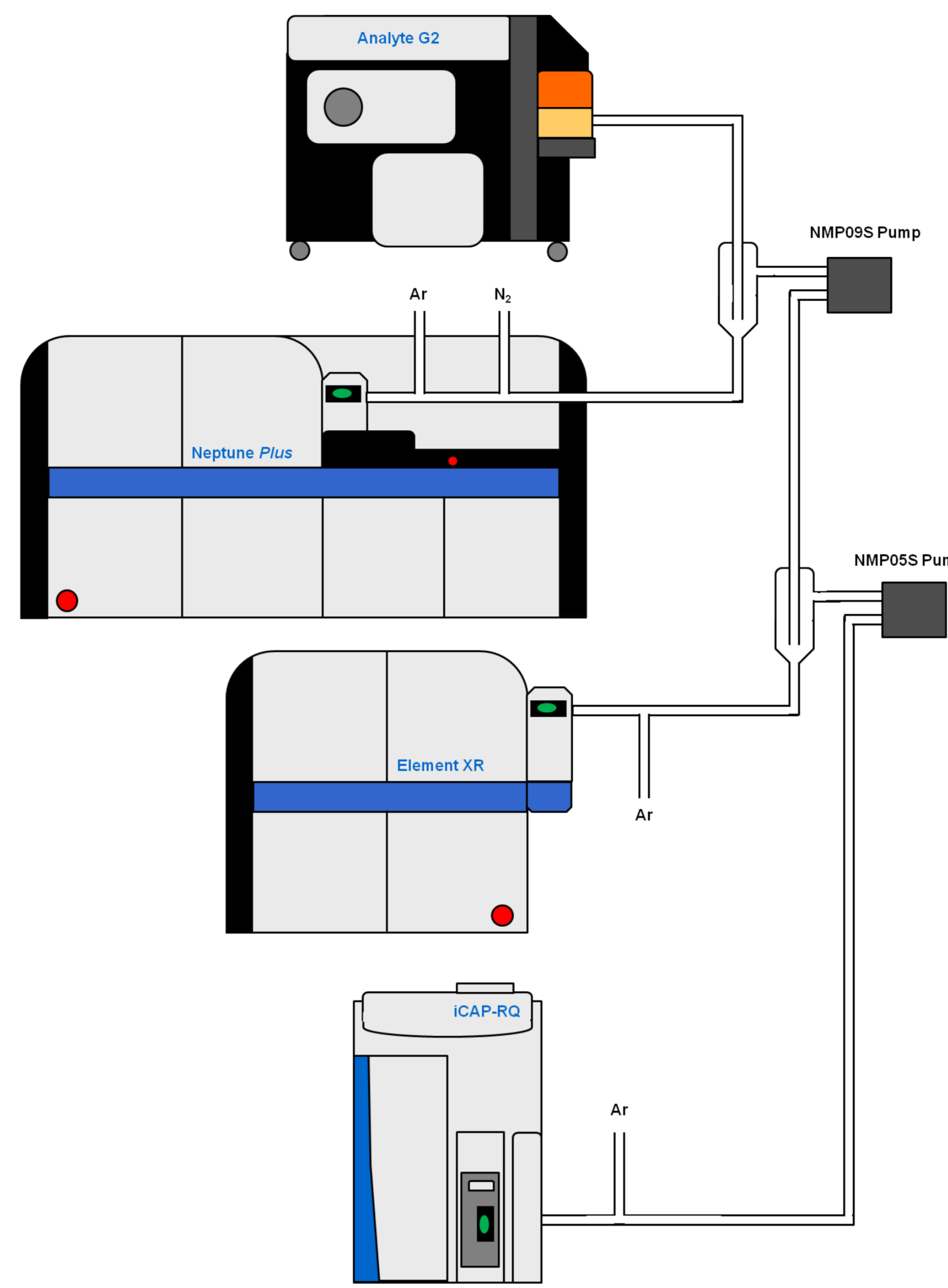


FIGURE 1. Schematic of LASS to three ICP-MS. Approx. 50% of the laser output was extracted from the primary MC-ICP-MS channel to SF-ICP-MS. Approx. 30% of the SF-ICP-MS channel extracted to Q-ICP-MS.

Results

Experimental

Each ICP-MS was tuned for high sensitivity, a U/Th ratio of approximately 1 and low oxides. The split was controlled via the pumps to supply sufficient signal on each ICP-MS during tuning.

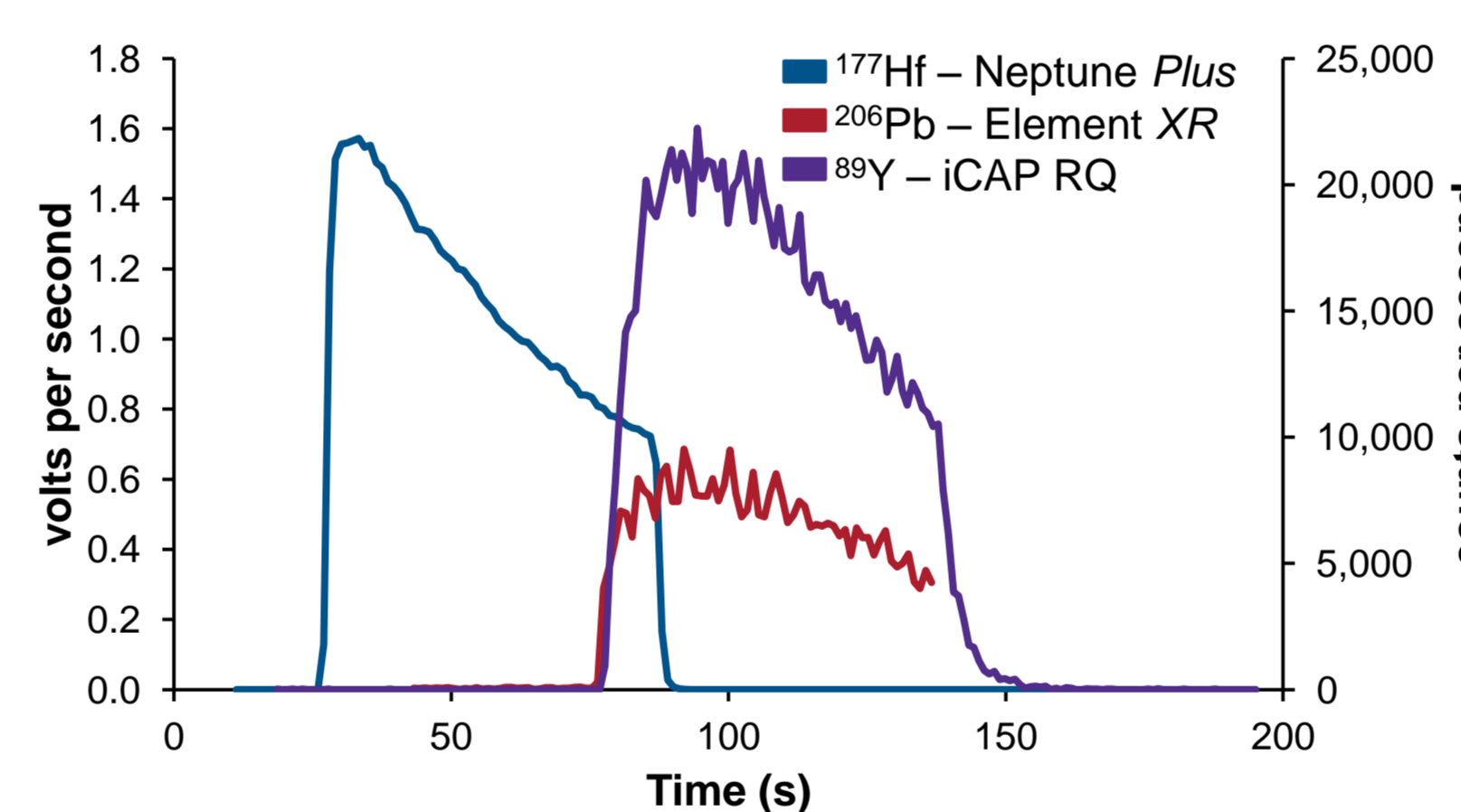


FIGURE 2. Signal traces for spot ablation of 91500 zircon, LASS to three ICP-MS. Integration time ≈ 1s. ¹⁷⁷Hf (MC-ICP-MS), responded faster than ²⁰⁶Pb (SF-ICP-MS) and ⁸⁹Y (Q-ICP-MS), reflecting the relative distances from the laser ablation cell.

Mean ¹⁷⁶Hf/¹⁷⁷Hf, U-Pb ages and REE concentrations were determined for 10 ablation craters. For REE analysis ⁹¹Zr was used as the internal standard. The 91500 zircon was used as an external standard, employing the sample-standard bracketing model. The other three reference materials, Plesovice, GJ-1 and Mud Tank, were used as unknown samples.

Plesovice

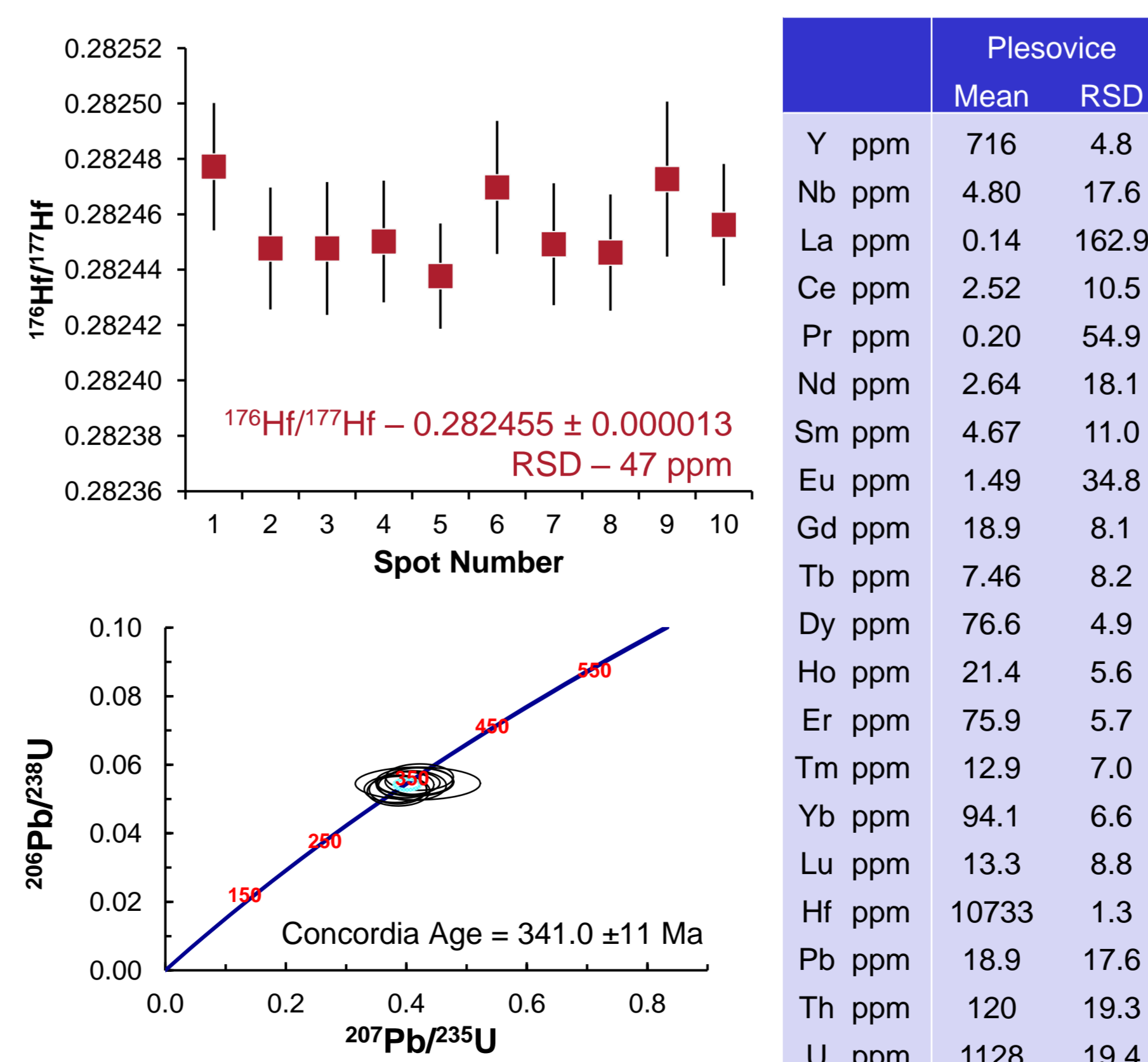


FIGURE 3. LASS results for Plesovice. ²⁰⁶Pb/²³⁸U Age – 339.9 ± 15.4 Ma (2σ).

GJ-1

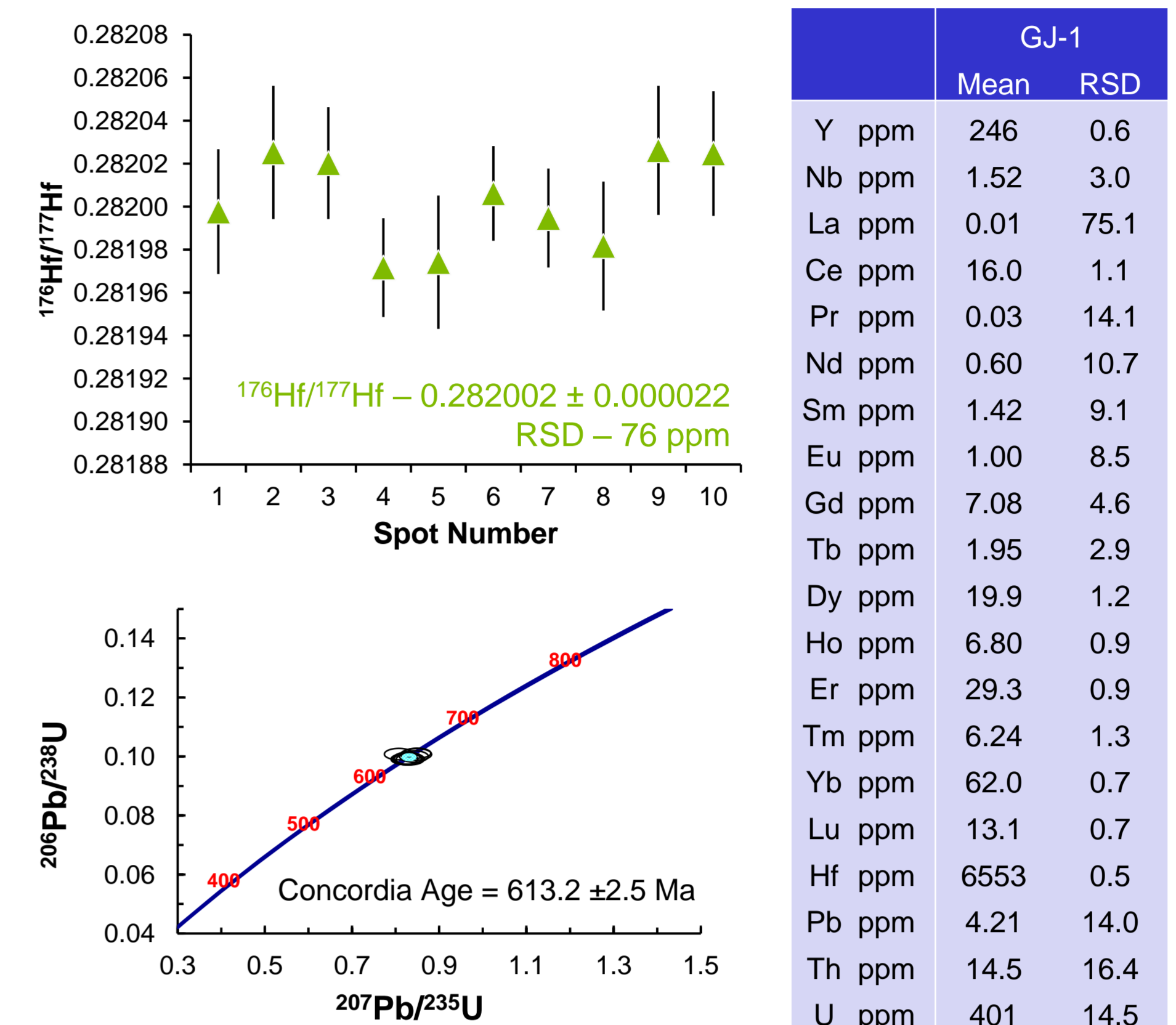


FIGURE 4. LASS results for GJ-1. ²⁰⁶Pb/²³⁸U Age – 612.5 ± 10.0 Ma (2σ).

Mud Tank

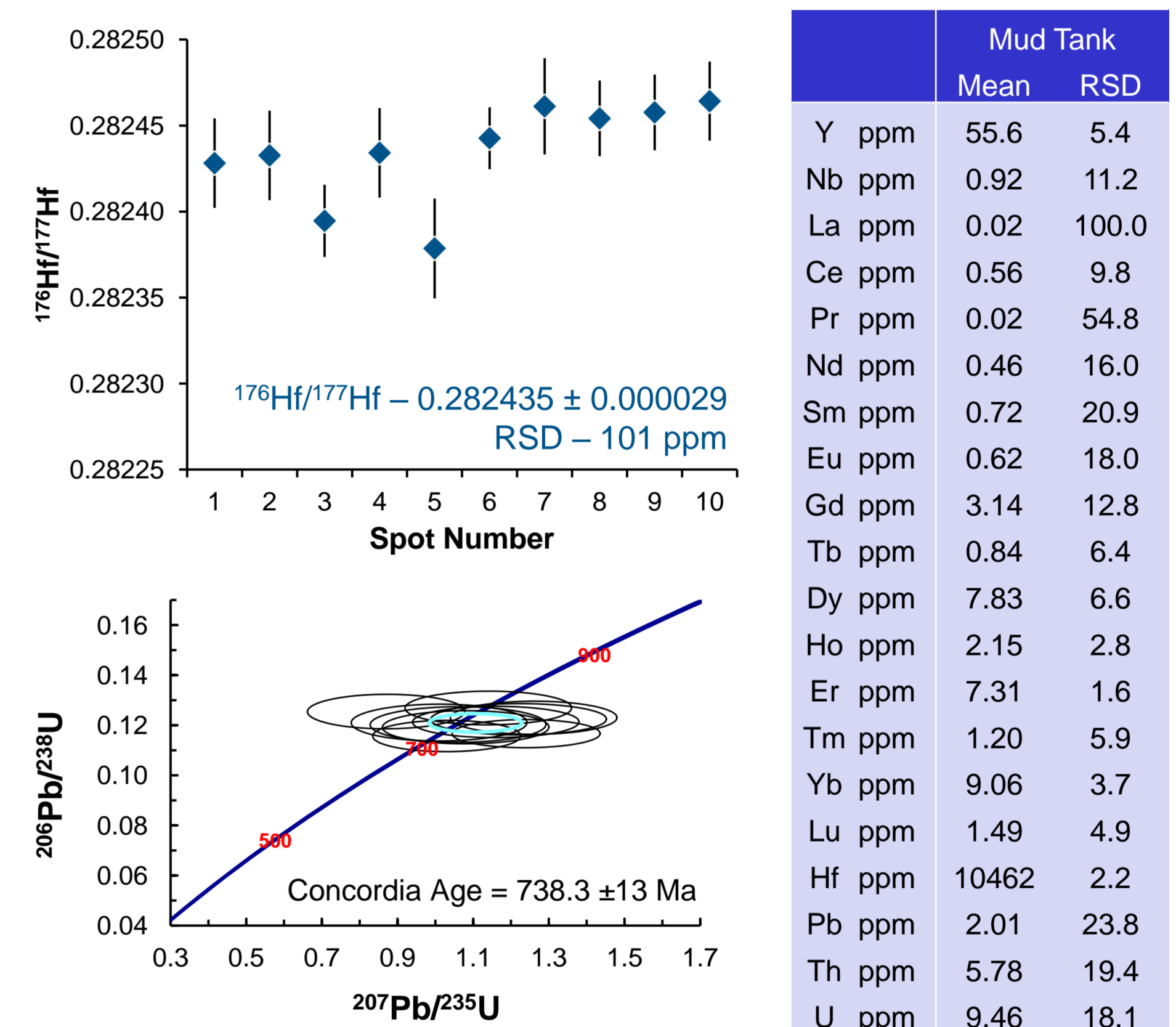


FIGURE 5. LASS results for Mud Tank. ²⁰⁶Pb/²³⁸U Age – 733.3 ± 39.9 Ma (2σ).

Hf

¹⁷⁶Hf/¹⁷⁷Hf ratios with RSD ≈ 100 ppm were achieved for all three sample reference materials.

U-Pb

The corresponding U-Pb ages were accurate, within uncertainty, to accepted values.

REE

The REE precision obtained varied depending on concentration and heterogeneity of the sample material.

Conclusion

The gas stream from a laser ablation system can be split between three ICP-MS.

- Membrane pumps can control the split to each ICP-MS
- Simultaneous Hf, U-Pb and REE determination

References

1. A. Kylander-Clark, B. Hacker, J. Cottle, *Chem. Geo.*, 2013, **345**, 99-112.
2. D. Frick, J. Schuessler, J. Buhk, F. von Blanckenburg, *Goldschmidt Abstracts*, 2015, **947**
3. C. Paton, J. Hellstrom, B. Paul, J. Woodhead, J. Hergt, *J. Anal. At. Spectrom.*, 2011, **26**, 2508-2518

Acknowledgements

"We acknowledge the IsoNose Initial Training Network funded by the People Programme (Marie Curie Actions) of the European Union's Seventh Framework Programme FP7/2007-2013/ under REA grant agreement number [608069]."

Teledyne Cetac Technologies, Teledyne Photon Machines, HelEx II and Analyte G2 are trademarks of Teledyne Instruments, Inc. Igor Pro is a trademark of Wavemetrics Inc. Iolite is a trademark of Iolite Software. SRM is a trademark of the National Institute of Science and Technology. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries.

This information is not intended to encourage use of these products in any manner that might infringe the intellectual property rights of others.

Thermo
SCIENTIFIC

A Thermo Fisher Scientific Brand